


File N:\Docs\Company\Manuals\D185\Accessories \ Output Connector.lwp  
Last Updated - March 30, 2004  
Printed - March 30, 2004

### Insertion of D185 Stimulus Output Connector

Please note that there are two levels of force needed to insert the Stimulus Output connector into the Stimulus Output Sockets on the D185.

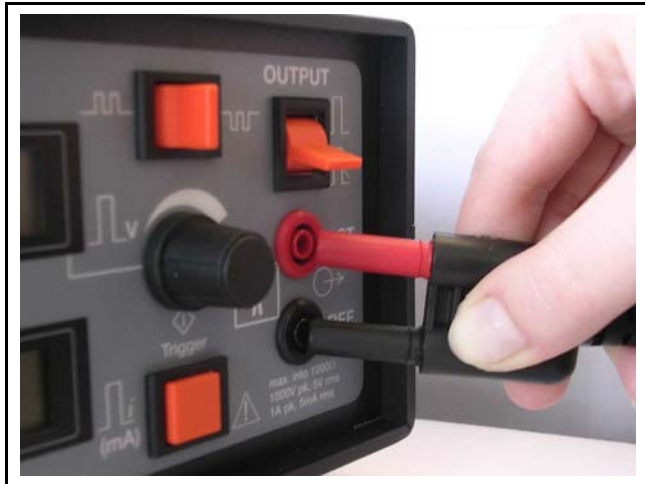

This photograph shows the correct orientation.

Red Plug to Red Socket.



To get to this point, very little force is required.

This photograph shows the connector **partially** inserted.

At this point increased resistance will be felt.



More force is required to fully insert the connector.

This zoomed photograph shows the connector **fully** inserted.