


D188 Remote Electrode Selector Stimulator Connection & Output Cable Options

The **D188 Remote Electrode Selector** is not supplied with stimulator connection cables or output extension cables. Input and output cabling requirements should be considered when ordering the D188.

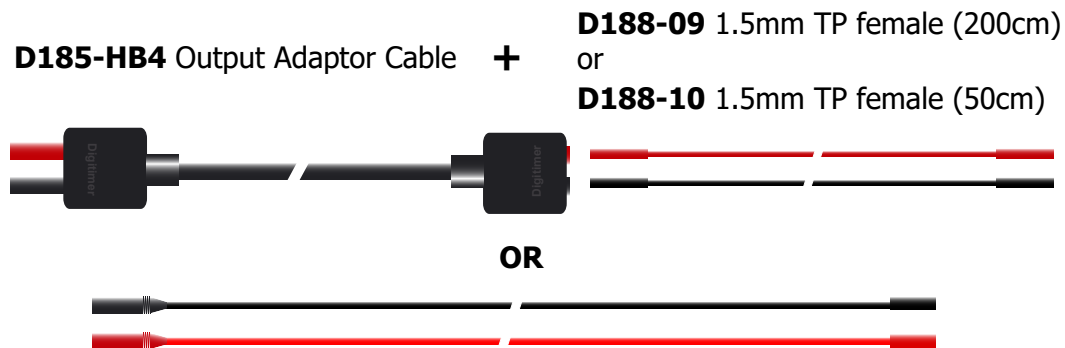

For direct connection from a Digitimer stimulator fitted with 4mm output sockets select **D188-12** 4mm TP male to 1.5mm TP female (50cm).

D188 INPUT

For connection to Digitimer stimulators with 2mm sockets (e.g. DS2A, DS3 or DS4), select the **D188-11** 2mm TP male to 1.5mm female (50cm).

D188 OUTPUT



Electrodes may be connected directly to the 1.5mm DIN42802 outputs on the D188.

In order to increase the distance between the subject and the D188 front panel or when electrodes leads are too short, pairs of 1.5mm TP D188 Output Extension Cables are available in 1m (**D188-13**) or 2m (**D188-14**) lengths.

Connection to D188 Output Sockets



Heatshrink is fitted at regular intervals to prevent cables from tangling.

Connection to electrode leads fitted with 1.5mm DIN42802 female connectors.